

Check Pesticide residues in fruits and vegetables with

Ringbio Rapid Test Kit










ltem#	Product
V001	Carbendazim Rapid Test Kit
V002	Paraquat Rapid Test Kit
V003	Chlorothalonil Rapid Test Kit
V004	Acetamiprid Rapid Test Kit
V005	Imidacloprid Rapid Test Kit
V006	Cyromazine Rapid Test Kit
V007	Carbaryl Rapid Test Kit
V008	Thiamethoxam Rapid Test Kit
V009	Triazolone-Triadimenol Test Kit
V010	Carbofuran Rapid Test Kit


ltem #	Product
V011	Albendazole Rapid Test Kit
V013	Fipronil Rapid Test Kit
V017	Carbendazim & Acetamiprid &
	Carbaryl 3in1 Rapid Test Kit
V018	Thiamethoxam & cyromazine, TC
	Combo Test Kit
V029	Glyphosate test kit
** 11.1.1	40 tt 00 tt b 0l- b#-

^{**} all kits are 48 tests or 96 tests per box. Sample buffer and test strips are included.


Procedures for testing

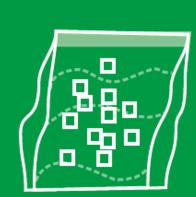

Take sample piece into a bag and mix with sample buffer

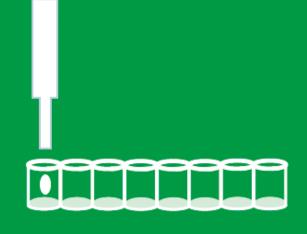

Add 200ul mixed sample into vial mix and incubate for 3min

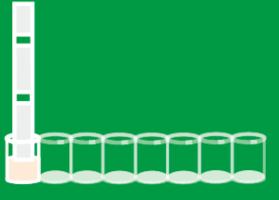

Insert test strip
Wait for 5min



Check the resu

^{**} all samples shall be homogenized before testing. For colored sample, filtration or centrifuge may require.

Please always check the kit instruction manual for detailed operation guide.